

# How to Install Patch 5

## ICP-MS MassHunter G7201A A.01.02


---

### Contents


|   |                                                                                                               |    |
|---|---------------------------------------------------------------------------------------------------------------|----|
| 1 | Uninstall G7201A A.01.01 (if it is installed).....                                                            | 2  |
| 2 | Install G7201A A.01.02 (if it is not installed). (skip this step and 3, if G7201A A.01.02 is installed) ..... | 6  |
| 3 | Updating System Parameters.....                                                                               | 7  |
| 4 | Install Patch 5 for ICP-MS MassHunter G7201A A.01.02.....                                                     | 8  |
| 5 | Confirm Build No. ....                                                                                        | 10 |

## How to Install Patch 5 for G7201A A.01.02

- 1 Uninstall G7201A A.01.01 (if it is installed)
  - 1.1 Click [Start] > [Control Panel] > [Programs and Features]

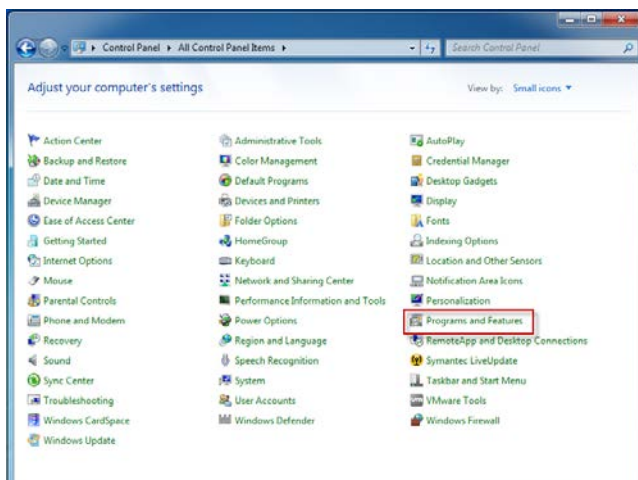


- 1.2 You will find "IcpLabQCAddin" and "ICP-MS MassHunter Workstation" in the table.

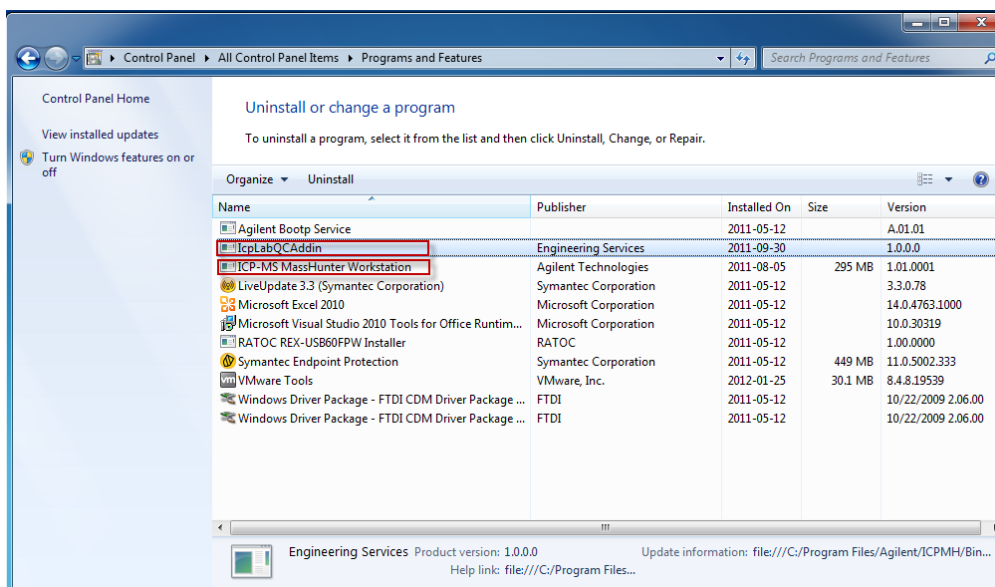


## How to Install Patch 5 for G7201A A.01.02

### 1.3 Right click on “IcpLabQCAddin” and select “Uninstall”

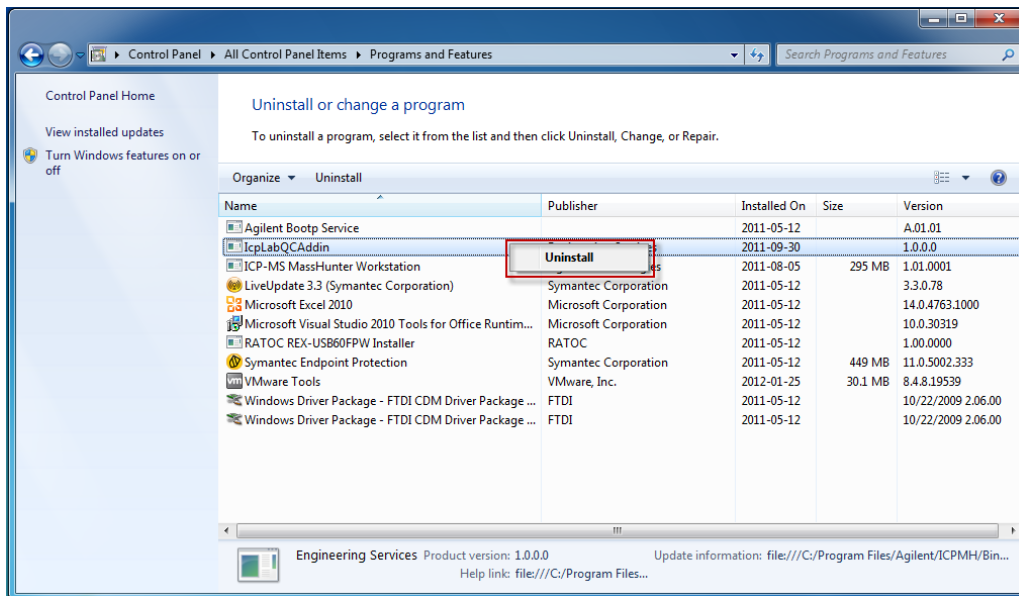


### 1.4 Right click on “ICP-MS MassHunter Workstation” and select “Uninstall”

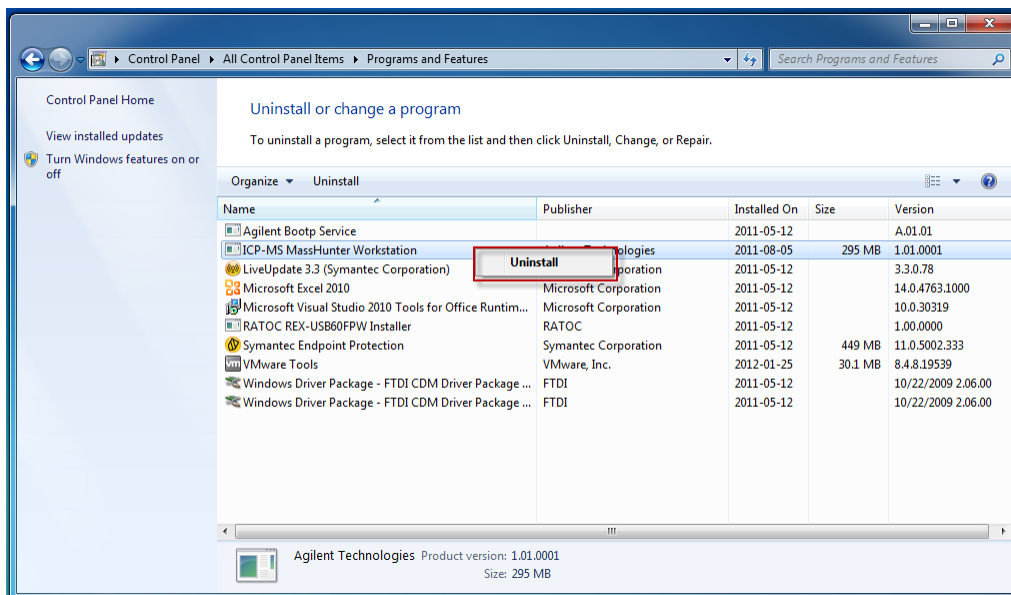


### 1.5 Close [Control Panel]

## How to Install Patch 5 for G7201A A.01.02

- 1.6 Open [Windows Explorer] and go to c:\Users\<user name>\AppData\Agilent\_Technologies,\_Inc

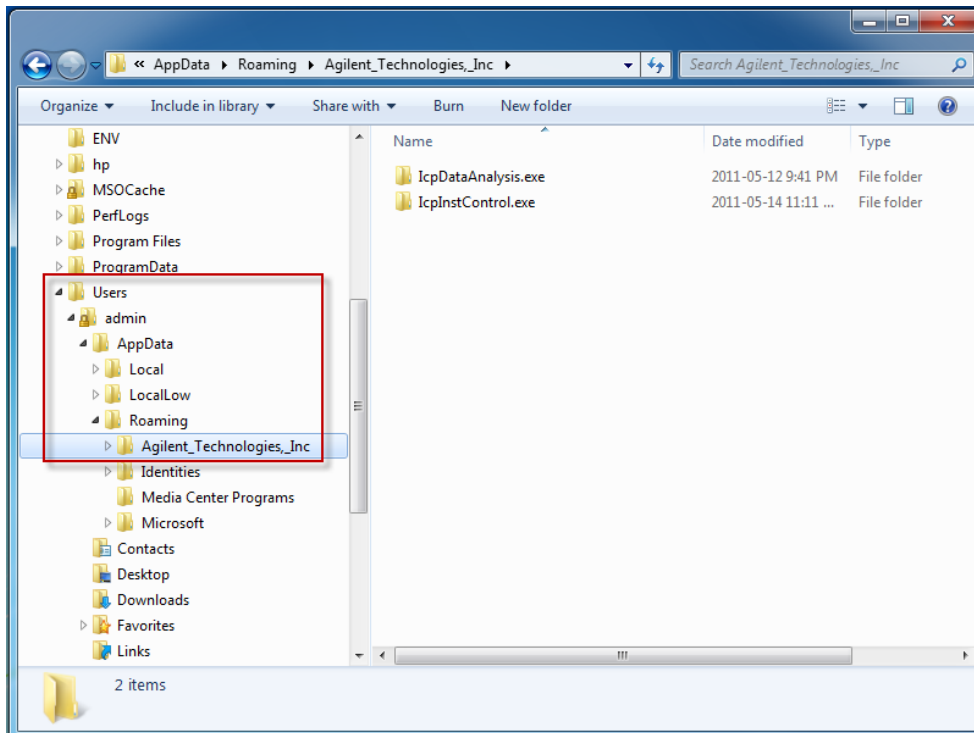


**Notice:** If you can't find the AppData folder, go to [Organize] > [Folder and Search options] > [View] tab > then select [Show hidden files, folders, and drives] under [Hidden files and folders].

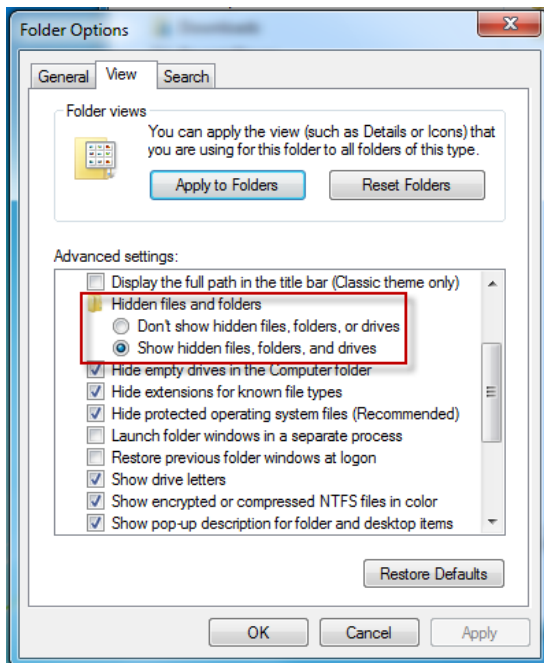


## How to Install Patch 5 for G7201A A.01.02

### 1.7 Delete the “IcpDataAnalysis.exe” and “IcpInstControl.exe” folders

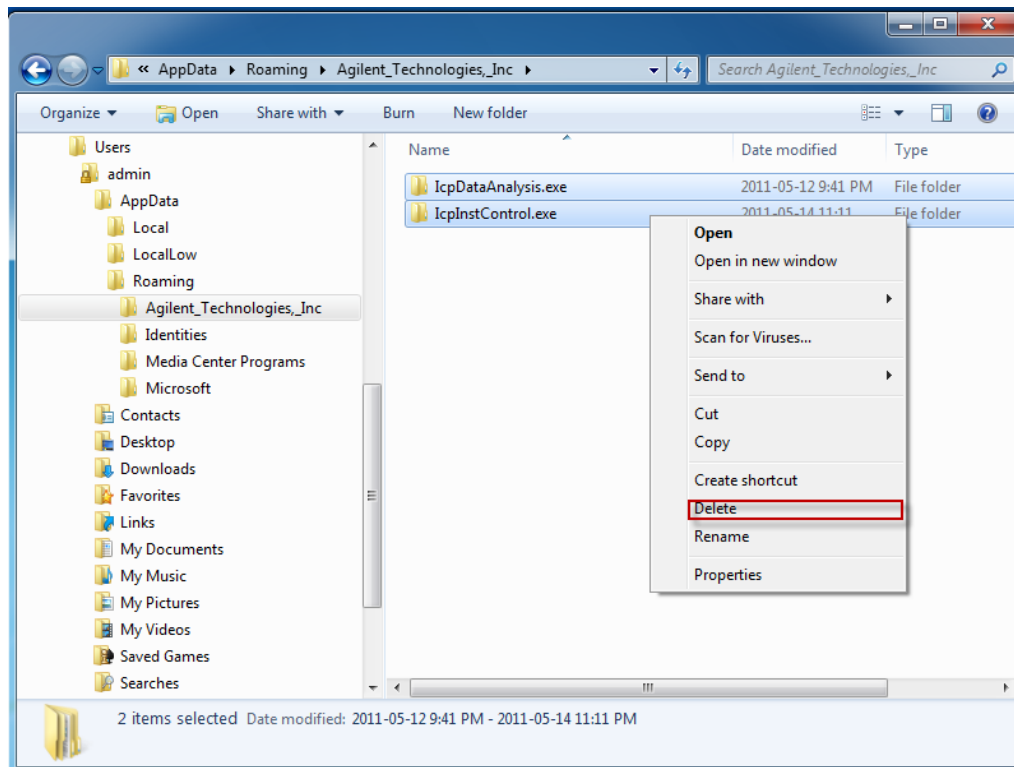


### 1.8 Delete “Configuration(.xml)” (in the path c:\agilent\ICPMH\1\)

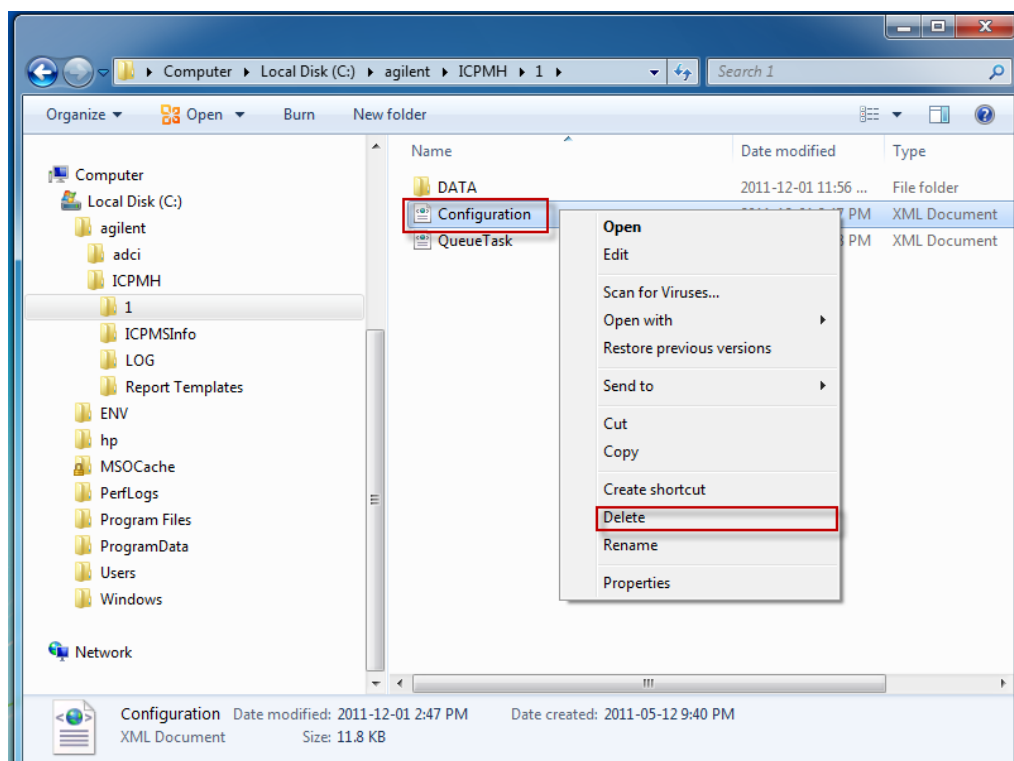

## How to Install Patch 5 for G7201A A.01.02

- 2 Install G7201A A.01.02 (if it is not installed). (skip this step and 3, if G7201A A.01.02 is installed)

- 2.1 Follow the instructions on the Index.htm page in the G7201A A.01.02 media DVD.

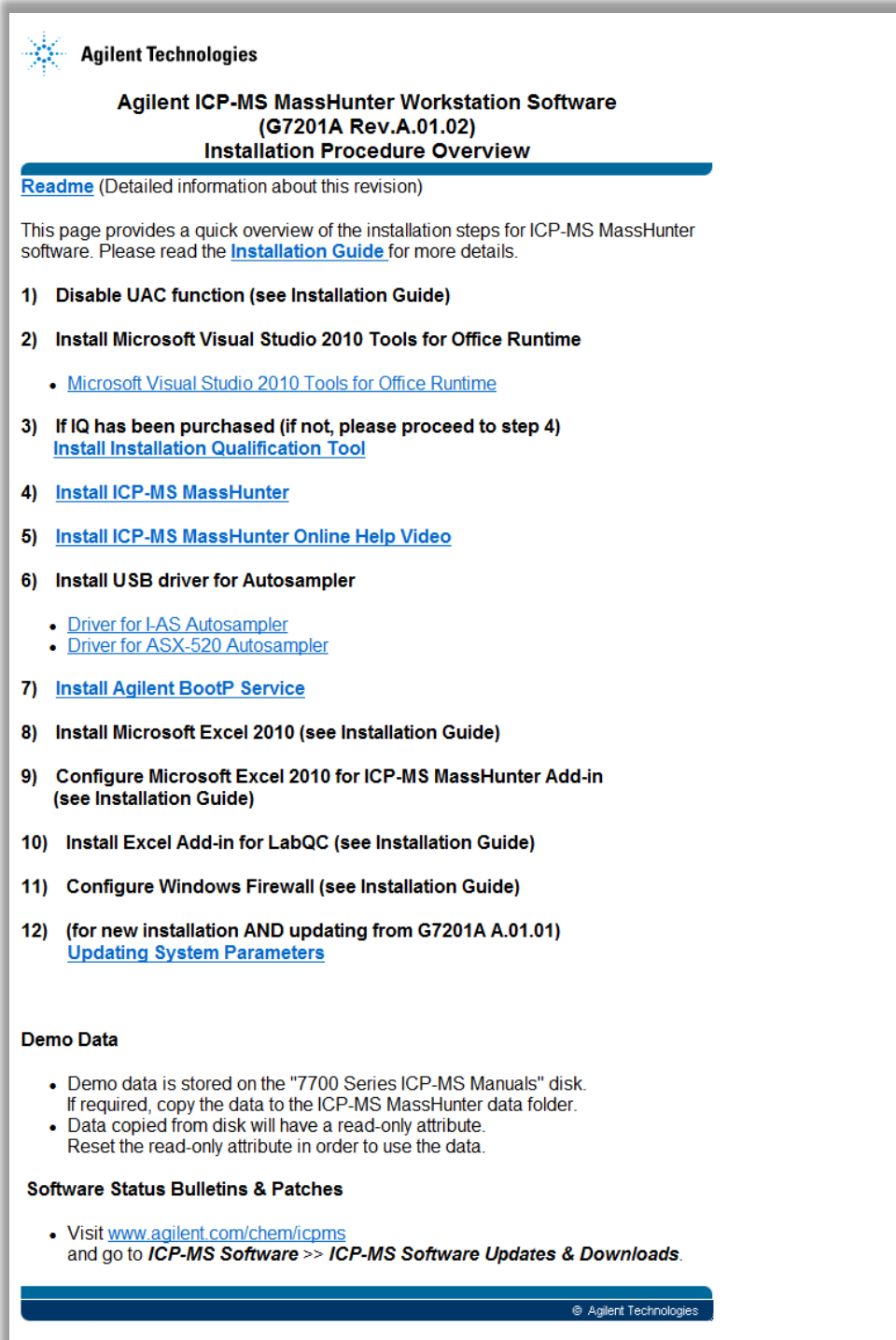

 **Agilent Technologies**

**Agilent ICP-MS MassHunter Workstation Software  
(G7201A Rev.A.01.02)  
Installation Procedure Overview**

[Readme](#) (Detailed information about this revision)

This page provides a quick overview of the installation steps for ICP-MS MassHunter software. Please read the [Installation Guide](#) for more details.

- 1) **Disable UAC function (see Installation Guide)**
- 2) **Install Microsoft Visual Studio 2010 Tools for Office Runtime**
  - [Microsoft Visual Studio 2010 Tools for Office Runtime](#)
- 3) **If IQ has been purchased (if not, please proceed to step 4)**  
[Install Installation Qualification Tool](#)
- 4) [Install ICP-MS MassHunter](#)
- 5) [Install ICP-MS MassHunter Online Help Video](#)
- 6) **Install USB driver for Autosampler**
  - [Driver for I-AS Autosampler](#)
  - [Driver for ASX-520 Autosampler](#)
- 7) [Install Agilent BootP Service](#)
- 8) **Install Microsoft Excel 2010 (see Installation Guide)**
- 9) **Configure Microsoft Excel 2010 for ICP-MS MassHunter Add-in (see Installation Guide)**
- 10) **Install Excel Add-in for LabQC (see Installation Guide)**
- 11) **Configure Windows Firewall (see Installation Guide)**
- 12) **(for new installation AND updating from G7201A A.01.01)**  
[Updating System Parameters](#)

**Demo Data**

- Demo data is stored on the "7700 Series ICP-MS Manuals" disk. If required, copy the data to the ICP-MS MassHunter data folder.
- Data copied from disk will have a read-only attribute. Reset the read-only attribute in order to use the data.

**Software Status Bulletins & Patches**

- Visit [www.agilent.com/chem/icpms](http://www.agilent.com/chem/icpms) and go to **ICP-MS Software >> ICP-MS Software Updates & Downloads**.

© Agilent Technologies

- 2.2 About 12) Updating System Parameters, refer to 3 Updating System Parameters.

### 3 Updating System Parameters


- 3.1 Some system parameters for the 7700 need to be updated by running the script (Ver2 Update System Parameter for ICP-MS MassHunterA0102.icpms.script).
- 3.2 Please refer to **Service Note 7700-031** (Plasma may not ignite with G7201A A.01.02) for instructions.
  - 3.2.1 Service Note  
[https://extranet.chem.agilent.com/sites/frontiersupport/\\_layouts/Agilent/HTMLVIEW/353089532.html](https://extranet.chem.agilent.com/sites/frontiersupport/_layouts/Agilent/HTMLVIEW/353089532.html)
  - 3.2.2 Instruction Guide  
[http://tad.jpn.agilent.com/support/Data/MassHunter\\_Workstation/Patches/Script/Instruction\\_for\\_updating\\_System\\_Parameters.pdf](http://tad.jpn.agilent.com/support/Data/MassHunter_Workstation/Patches/Script/Instruction_for_updating_System_Parameters.pdf)
- 3.3 The script can be found at either of following sites
  - 3.3.1 [http://tad.jpn.agilent.com/support/Data/MassHunter\\_Workstation/Patches/Script/New%20Script%20\(Ver2\).zipx](http://tad.jpn.agilent.com/support/Data/MassHunter_Workstation/Patches/Script/New%20Script%20(Ver2).zipx)
  - 3.3.2 MassHunter Workstation Software DVD (G7201A Revision A.01.02 **February 2012, G7201-60502**)

## How to Install Patch 5 for G7201A A.01.02


### 4 Install Patch 5 for ICP-MS MassHunter G7201A A.01.02

**Notice: Patch 5 includes Patch 001, Patch 002, Patch 003, and Patch 004 contents. It is not necessary to install Patch 001, Patch 002, Patch 003, or Patch 004 when installing Patch 5. Patch.msp includes all fixes for English, Chinese, and Japanese versions.**

- 4.1 Put the 7700 in shutdown mode from MassHunter.
- 4.2 Close MassHunter.
- 4.3 Copy patch file **“Patch.msp”** to the MassHunter PC.
- 4.4 Double click the patch file.
- 4.5 Click **[Next]**.

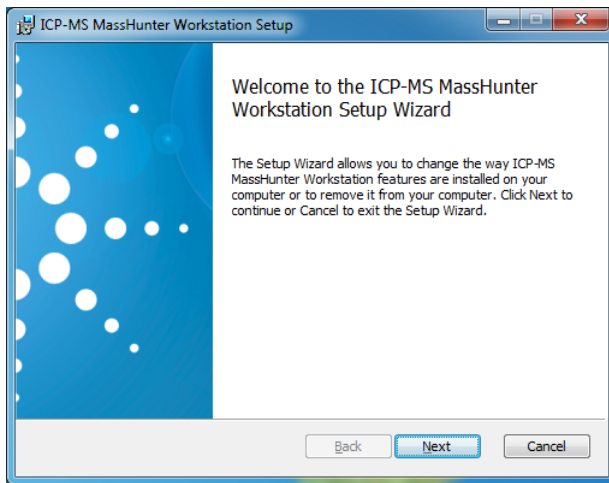


- 4.6 Click **[Repair]** and **[Next]**.

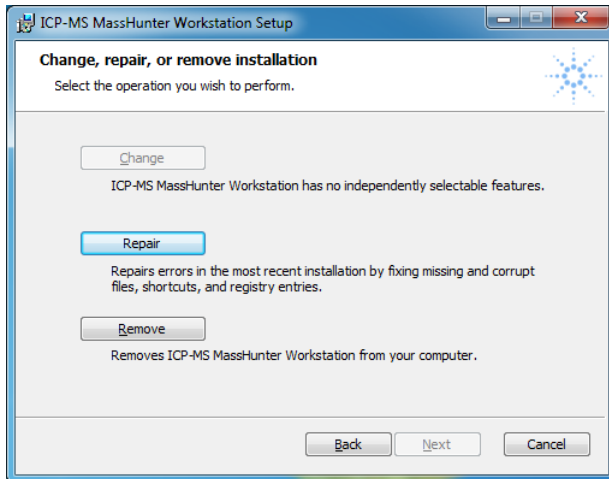


## How to Install Patch 5 for G7201A A.01.02

### 4.7 Click **[Repair]**.

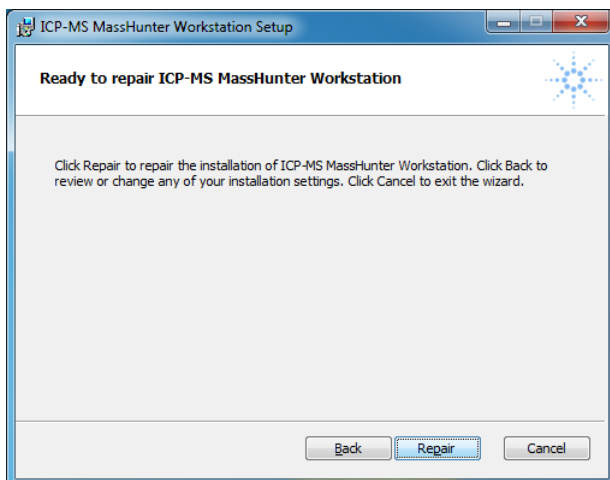


### 4.8 Click **[Finish]**.

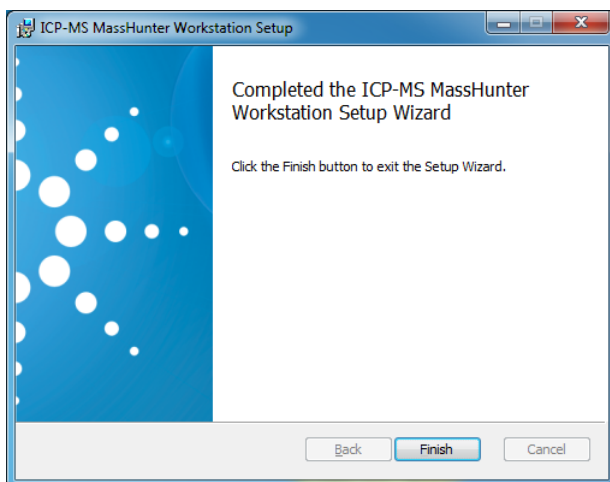

## How to Install Patch 5 for G7201A A.01.02

### 5 Confirm Build No.

- 5.1 Open **[Offline Data Analysis]**.
- 5.2 Select **[Help] > [About...]**, and the splash screen will appear indicating the Version, Build number and Patch number as follows (Version A.01.02 G7201A Build 291.22 Patch 5).

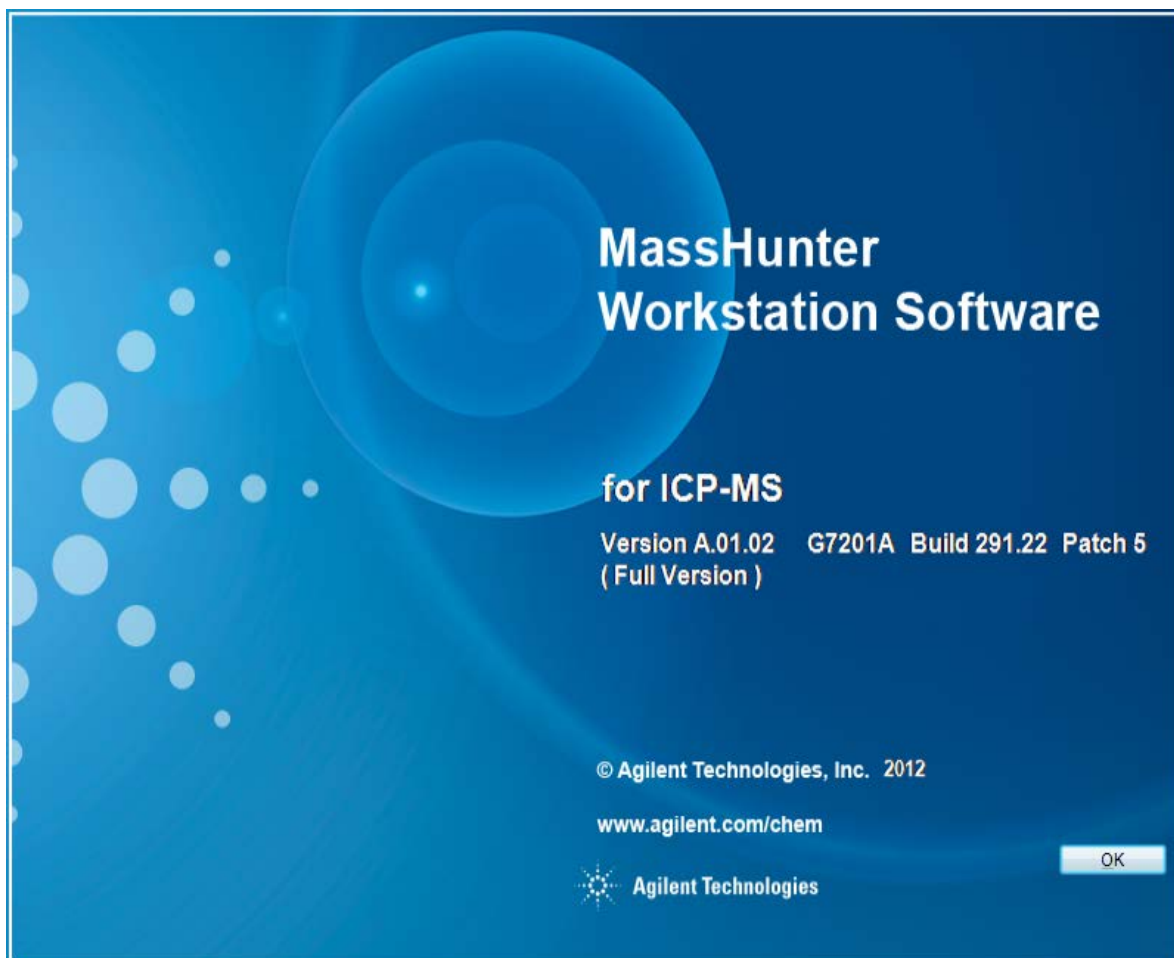

This information is subject to change without notice

© Agilent Technologies, Inc. 2012

Published in Japan, 01-October-2012