



Page 1 of 2

SERVICE How To

Model Number:	Originator:	Topic
PS 335 DAD	Des Wichems	Typing in the Optics Calibration Coefficients into the Firmware

PS 335 Transferring the Calibration Coefficients in to the Firmware

- [illegible]

2. Open Diagnostics and go to the *Calibration* tab (1) (see screen capture below).
3. You can display the current calibration settings by selecting the *Get Calibration Data* (2) option. Once they are displayed you will need to select the *Write* (3) option and then type in the wavelength coefficients (on the sticker on the top of the new optics module) into the shown calibration file.
4. Once this is completed, you will need to save (4) the new coefficients, and then you can then do a lamp cal in either Star or Galaxie software.

