D.03.00 SP2 *MUST* be installed onto a previously installed D.03.00 SP1.

If you have installed any point patches (e.g. .dll or .mac files) onto your D.03.00 SP1, that is OK; you do not have to remove them or do anything special before you install D.03.00 SP2.

After installation, consult the file X:\msdchem\msexe\ReadmeD0300sp2.txt for a list of the defects fixed in this revision.

Important Design Changes:

The software repair for PVCS #1076 required a design change to the Sample Log Table (Sequence Editor). The previous design would allow methods and data file names to be selected from a file browser. The file browser would be invoked when the question mark (“?”) was selected in the Method or Data File column. The current (new) design invokes the same file browser, but from a right-click pop-up menu item on a cell in the Method or Data File column.

Also, when you are using a Keyword “MethPath” or “DataPath”, you no longer have to manually enter method or data file paths into the KeywordString column. Now, you can right-click in the KeywordString column and use a file browser to select your method or data file paths.

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