

$$\Theta = \pm 45^\circ \frac{\lambda^2}{\lambda_0^2} \left(\frac{\lambda_0^2 - F}{\lambda^2 - F} \right)^2$$



Laser Modulation



$$\Theta = V \int_{z_1}^{z_1+L} H_1 dz$$



Introduction

Conoptics manufactures an extensive line of low voltage electro-optic light modulators, driver electronics, and associated components to satisfy diverse requirements. The application dictates which versions of modulator and driver electronics are to be used and what auxiliary components are needed. This data sheet presents the characteristics of standard products and their operating parameters when combined to form functional systems. Laser noise reduction systems, versions of modulation systems with special feedback electronics, are described in separate data sheets. In addition to catalog products, Conoptics also manufactures, on a custom basis, intensity modulation systems with high frequency capabilities beyond 400 MHz and frequency chirping and FM spectroscopy modulators out to several GHz.

Drive Electronics

In general, the first application requirements considered in the choice of modulation system components are the information bandwidth and waveform requirement. The driver output voltage achievable is a function of amplifier bandwidth and, while this system parameter is not isolated from others, such as aperture diameter, operating wavelength, etc., it is normally the limiting parameter of the system. Standard Conoptics products include four general purpose drivers:

The Models 10, 25, 50 and 100 are dc coupled broadband amplifiers which require an input of 1 volt p-p into 50 ohms for full rated output. Their bandwidths are a function of the modulator used. Each model requires a different electrical configuration in the associated modulator. The Model 10 requires that the modulator be configured as a lumped capacitor. The Model 25 requires a 100 ohm balanced line; the Model 50, a 50 ohm balanced line;; and the Model 100, a two segment (4 port) 50 ohm balanced line.

POLYTEC GmbH

Büro Berlin Schwarzschildstraße 1 D - 12489 Berlin GERMANY
Tel: +49 (30) 63 92 51 40 Fax: +49 (30) 63 92 51 41



wl@polytec.de
www.polytec.de

$$\Theta = \pm 45^\circ \frac{\lambda_z^2}{\lambda_o^2} \left(\frac{\lambda_o^2 - F}{\lambda^2 - F} \right)^2$$



Laser Modulation



$$\Theta = V \int_z^{z+L} H_x dz$$

The Model 302 is primarily intended for low signal bandwidth, long optical wavelength applications. It also offers cost advantages over higher frequency broadband drivers, especially since, due to its high voltage output, shorter capacitive modulators can be used. The bandwidth of the Model 302 ranges from 150 to 200 KHz depending on the modulator used. Input requirement is 4 volts p-p into 50 ohms. All models include a built in manual bias control.

Optical Modulators

All modulators listed in this data sheet are of the transverse field type, that is, the electric field produced by the applied signal voltage is perpendicular to the optical propagation direction. The voltage swing required by a given modulator at a given operating wavelength to transit between the full off state to the full on state is called the Half Wave Voltage ($V_{\frac{1}{2}}$). The transverse field structure allows reduction of $V_{\frac{1}{2}}$ by manipulation of the crystal length to aperture ratio to a level achievable by available driver electronics. $V_{\frac{1}{2}}$ is roughly proportional to wavelength and long wavelength devices usually require higher length to aperture ratios to accommodate existing driver output levels.

Conoptics offers modulators constructed with three different crystal species: Ammonium Dihydrogen Phosphate (ADP), Potassium Dideuterium Phosphate (KD*P), and Lithium Tantalate (LTA).

Models 370, 380, and 390 utilize ADP as the active element. The unique feature of these models is the virtual non-existence of piezoelectric resonances.

Models belonging to the 360 Series utilize LTA. LTA has the lowest intrinsic $V_{\frac{1}{2}}$ and the longest wavelength IR cutoff. Furthermore, it has a combination of high refractive index and relatively low dielectric constant which allows modulators to be designed which make full use of the intrinsic driver frequency response. Models in the 360 Series exhibit piezoelectric resonances but they are discrete and very narrow.

KD*P is used in Model 350 Series modulators. In terms of optical transmission bandwidth and driver frequency response utilization, this series falls in between ADP and LTA versions.

$$\Theta = \pm 45^\circ \frac{\lambda^2}{\lambda_0^2} \left(\frac{\lambda_0^2 - F}{\lambda^2 - F} \right)^2$$



Laser Modulation



$$\Theta = V \int_z^{z+L} H_1 dz$$

Modulator Modifications

Any of the modulators listed here can be used as a phase modulator by simply rotating the input polarization direction by 45°. This procedure makes one of the modulator half segments essentially inactive and doubles $V^{1/2}$ (now the voltage required for a 180° phase shift). A factory modification can be made during construction which restores $V^{1/2}$ to its original value. This modification precludes use of the device as an intensity modulator, however, and is irreversible.

Auxiliary Components

With the exception of 360 Series, modulators used at wavelengths longer than 2000nm, an integral Glan type polarizer (analyzer) is supplied with each model listed here. Operation at longer wavelengths requires polarizers of a different type and may be additional cost items. Other components such as quarter wave plates used in polarization rotators, are also available from Conoptics.

The most commonly used auxiliary components are Automatic Bias Controllers (ABC's). The purpose of ABC's is to compensate the long term temperature induced drift of the bias voltage needed to position the applied signal baseline at the desired operating point on the modulator transfer characteristic. Three different versions are available. The first accommodates signal information flows which have a periodic "dead time" such as scanned data or that found in image recorders. Here, a sampling signal is injected by the ABC during the "dead time" and the resulting optical modulation is analyzed to produce an error signal. A feedback loop drives the operating point to the top or bottom of the transfer characteristic, as desired. The second option, used with continuous information flows, such as video disc mastering, samples both the modulated optical output and its reciprocal signal. It averages these samples and produces an error signal which drives the operating point to the midpoint of the transfer characteristic. The third option is similar to the second but is designed to control arbitrary duty cycle digital waveforms.

All ABC versions are available with modulation systems incorporating ADP or KD*P modulators and Model 10, 25, 50 and 100 driver electronics. The inherent stability of 360 Series LTA modulators is sufficient in the majority of applications to avoid the need for an ABC. The addition of an ABC to a modulation system requires integration with both the driver electronics and the optical modulator and is a factory installed option.

POLYTEC GmbH

Büro Berlin Schwarzschildstraße 1 D - 12489 Berlin GERMANY
Tel: +49 (30) 63 92 51 40 Fax: +49 (30) 63 92 51 41



wl@polytec.de
www.polytec.de

$$\Theta = \pm 45^\circ \frac{\lambda^2}{\lambda_0^2} \left(\frac{\lambda_0^2 - F}{\lambda^2 - F} \right)^2$$



Laser Modulation



$$\Theta = \int_{z_0}^{z_0+L} H_1 dz$$

Modulation Systems

The modulators and drivers listed in this data sheet can be used in various combinations to form high performance, cost effective modulation systems.

Table II shows the key performance characteristics of various combinations of standard driver electronics and modulators. The high frequency -3dB points may be limited either by the driver or the modulator. Rise and fall times are normally calculated as 0.35 divided by the -3dB bandwidth but, due to the compression caused by the sine squared transfer characteristic over its full on to off range, the optical rise and fall times of these systems is approximately 20% less.

Table 1 Modulator Specifications:

Note: All cells are 50mm diameter.

Model No	350-50	350-80	350-105	350-160	350-210	360-40	360-80	360-120	360-160	370	370 LA	380	390
Crystal	KD*P	KD*P	KD*P	KD*P	KD*P	LTA	LTA	LTA	LTA	ADP	ADP	ADP	ADP
V 1/2 , volts @ 500 nm 830 nm 1064nm 2500nm	455 757 NA NA	261 433 522 NA	226 376 482 NA	130 216 NA NA	113 188 NA NA	NA 312 400 NA	NA 143 183 430	NA 107 138 323	NA 71 92 215	184 306 NA NA	263 437 NA NA	92 153 NA NA	115 190 NA NA
Aperture Diameter, mm	3.1	2.7	3.1	2.7	3.1	2.7	2.7	2.7	2.7	2.5	3.5	2.5	3.5
Useful Transmission Range, nm	300- 1100	300- 1100	300- 1100	300- 1000	300- 1000	700- 4500	700- 4500	700- 4500	800- 4500	300- 750	300- 750	300- 750	300- 750
Resonances	YES	YES	YES	YES	YES	YES	YES	YES	YES	NO	NO	NO	NO
Contrast Ratio @ 633nm 1064nm	500:1 700:1	500:1 700:1	500:1 700:1	300:1 NA	300:1 NA	NA 300:1	NA 200:1	NA 200:1	NA 100:1	500:1 NA	500:1 NA	500:1 NA	500:1 NA
Length, mm with Polarizer	106	137	162	215	268	95	137	174	215	158	158	253	272

POLYTEC GmbH

Büro Berlin Schwarzschildstraße 1 D - 12489 Berlin GERMANY
Tel: +49 (30) 63 92 51 40 Fax: +49 (30) 63 92 51 41



wl@polytec.de
www.polytec.de

$$\Theta = \pm 45^\circ \frac{\lambda^2}{\lambda_0^2} \left(\frac{\lambda_0^2 - F}{\lambda^2 - F} \right)^2$$



Laser Modulation



$$\Theta = \int_{z_0}^{z_0+L} H_1 dz$$

Table 2 Modulation Systems:

MODULATOR MODEL NUMBER	10	25	50	100	200	302	302A	25D
350-50	*	*	*	*	1 to 200 MHz	DC - 200KHz 875nm	DC-1MHz	*
350-80	*	*	*	*	1 to 200 MHz	CD - 200KHz 1100nm	DC-1MHz	<8ns,>25MHz 354nm
350-105	*	*	*	*	1 to 140MHz	*	DC-1MHz	<8ns,>20MHz 430nm
350-160	DC - 12MHz 460nm	DC - 35MHz 460nm	*	*	1 to 100 MHz	*	DC-1MHz	<8ns,>20MHz 710nm
350-210	DC - 12MHz 530nm	DC - 30MHz 530nm	DC - 60MHz 400nm	CD - 120MHz 400nm	1 to 80MHz	*	DC-1MHz	<8ns,>20MHz 820nm
360-40	*	*	*	*	1 to 200 MHz	DC - 300KHz 2120nm	DC-1MHz	*
360-80	*	*	*	*	1 to 200 MHz	DC - 300KHz 4500nm	DC-1MHz	<8ns,>20MHz 1010nm
360-160	DC - 13MHz 1390nm	DC - 40MHz 1390nm	DC - 80MHz 1040nm	DC - 130MHz 1040nm	1 to 125 MHz	*	DC-1MHz	<8ns,>20MHz 2135nm
360-120	DC - 12MHz 2090nm	DC - 35MHz 2090nm	DC - 60MHz 1575nm	DC - 120MHz 1575nm	1 to 150 MHz	*	DC-1MHz	<8ns,>10MHz 3235nm
370	*	*	*	*	1 to 125 MHz	DC - 200KHz 850nm	DC-1MHz	<8ns,>20MHz 505nm
370 LA	*	*	*	*	1 to 125 MHz	DC - 200KHz 850nm	DC-1MHz	<8ns,>20MHz 355nm
380	DC - 10MHz 650nm	DC - 25MHz 650nm	DC - 50MHz 490nm	DC - 100MHz 490nm	1 to 80 MHz	*	DC-1MHz	<8ns,>20MHz 80nm (2)
390	DC - 10MHz 520nm	DC - 25MHz 520nm	DC - 50MHz 400nm	DC - 100MHz 400nm	1 to 80 MHz	*	DC-1MHz	<8ns,>10MHz 810nm

* Configuration not typical. Call factory for depth of modulation.

POLYTEC GmbH

Büro Berlin Schwarzschildstraße 1 D - 12489 Berlin GERMANY
Tel: +49 (30) 63 92 51 40 Fax: +49 (30) 63 92 51 41



wl@polytec.de
www.polytec.de